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## SPEX CertiPrep®

## Certificate of Analysis

As you know, every manufacturer of Certified Reference Materials supplies a certificate of analysis with their product. Not all certificates are alike. We know because SPEX CertiPrep® has been supplying the most comprehensive certificate in the industry for years. Many other companies have followed, but no one gives you the information you get from us. We have highlighted why our certificate is the best and what you should look for in a certificate of analysis.

SPEX CertiPrep is accredited by A2LA for Organic and Inorganic Certified Reference Materials as complying with the requirements of ISO/IEC 17025 and ISO/IEC Guide 34 with the most comprehensive scope in the industry.

**SPEX Certificate**  
Certificate of Reference Material  
**CLARITAS PPT**

Lot No. CL6-154YP

**Catalog Number:** CL-ICS-1  
**Description:** Instrument Check Standard 1  
**Matrix:** 2% HNO<sub>3</sub> / Tr. Tart. Acid / Tr. HF (Ultra Pure)

This CLARITAS PPT® Certified Reference Material, CRM, is intended primarily for use as a calibration standard or quality control standard for inorganic spectroscopic instrumentation such as ICP-OES, DCP, AA, ICP-MS, and XRF. It can be employed in USEPA, ASTM and other methods relevant to the certified properties listed below. The CRM is prepared from high purity single element concentrates of individual elements using Class A laboratory ware to give precise concentrations. See side 2 for details of certification.

**Instrumental Analysis by ICP Spectrometer:**

Analyte	Labeled	Certified	Uncertainty	SRM	Analyte	Labeled	Certified	Uncertainty	SRM
Ag	10 µg/mL	9.90 µg/mL	±0.05 µg/mL	3151*	Mn	10 µg/mL	9.85 µg/mL	±0.05 µg/mL	3152*
Al	10 µg/mL	9.88 µg/mL	±0.05 µg/mL	3101a*	Ni	10 µg/mL	9.98 µg/mL	±0.05 µg/mL	3136*
As	10 µg/mL	9.91 µg/mL	±0.05 µg/mL	3103a*	Pb	10 µg/mL	9.95 µg/mL	±0.05 µg/mL	3130*
Ba	10 µg/mL	9.93 µg/mL	±0.05 µg/mL	3104a*	Sb	10 µg/mL	9.83 µg/mL	±0.05 µg/mL	3102a*
Be	10 µg/mL	9.87 µg/mL	±0.05 µg/mL	3105a*	Se	10 µg/mL	9.94 µg/mL	±0.05 µg/mL	3149*
Cd	10 µg/mL	9.86 µg/mL	±0.05 µg/mL	3108*	Tl	10 µg/mL	9.87 µg/mL	±0.05 µg/mL	3158*
Co	10 µg/mL	9.94 µg/mL	±0.05 µg/mL	3113*	V	10 µg/mL	9.87 µg/mL	±0.05 µg/mL	3185*
Cr	10 µg/mL	9.81 µg/mL	±0.05 µg/mL	3112a*	Zn	10 µg/mL	9.94 µg/mL	±0.05 µg/mL	3188a*
Cu	10 µg/mL	9.82 µg/mL	±0.05 µg/mL	3114*					

\* - indicates NIST SRM  
\* - indicates SPEX CertiPrep CRM (when NIST SRM is not available)

SPEX CertiPrep Reference Multi: Lot# 26-19JB-6-293GS-42-32AS

**Density:** 1.010 g/mL @ 20°C

**Trace Metallic Impurities in the Actual Solution via ICP-MS Analysis:**

Element	µg/L	Element	µg/L	Element	µg/L	Element	µg/L				
Au	<10	Fe	<3	K	<300	P	<700	Sc	<0.1	Th	0.2
B	1	Ga	0.8	La	1	Pd	<0.3	Si	<100	Ti	<0.7
Bi	0.5	Gd	<0.01	Li	<0.3	Pr	0.05	Sm	0.09	Tm	<0.01
Ca	<10	Hf	<0.3	Lu	<0.01	Pt	<0.06	Sr	<0.1	U	0.05
Ce	0.2	Hg	0.3	Mg	0.9	Rb	<0.09	Ta	2	W	<0.8
Cs	<0.1	Ho	<1	Mo	<0.5	Rh	<0.04	Tb	2	Y	0.03
Dy	<0.02	In	<0.01	Nb	0.5	Ru	0.5	Tc	<0.01	Yb	<0.01
Er	<0.02	Ir	<0.05	Nd	<0.02					Zr	0.7
Eu	0.8	Y	<0.03								

Date of Certification: \_\_\_\_\_  
Certifying Officer: \_\_\_\_\_  
Quality Officer: \_\_\_\_\_

68 elements are scanned with *found values* for Claritas PPT® and Assurance® Standards.

Each elemental impurity listed with *actual value* - not limited to elements above detection limits.

Trace impurities of the *final solution* - not of the starting material.

Stability and accuracy of the *final solution* - not the starting materials.

Traceable to NIST

Signed by SPEX CertiPrep®'s Vice President of Manufacturing and Quality Officer.

Stamped with month and year of shipment.

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## Report of Certification

This Certified Reference Material (CRM) has been prepared and certified under an ISO 9001:2000, ISO 17025:2005, and ISO Guide 34:2000 quality system consistent with the following guides:

- Guide To The Expression Of Uncertainty In Measurement 1997
- EURACHEM/CITAC Guide: Quantifying Uncertainty in Analytical Measurement – Second Edition
- ASTM Guide D6362-98
- ISO Guide 31: Contents of certificates of reference materials
- ISO Guide 34: Quality system guidelines for the production of reference materials – certified by A2LA
- ISO 17025: Certification of reference materials, general and statistical principles – certified by A2LA
- NIST Technical Note 1297
- ILAC-G12-2000: Guidelines for the requirements for the competence of reference materials producers
- ISO/REMCO N280

### Material Source:

All analytes and matrix materials are obtained and verified by SPEX CertiPrep from pre-qualified vendors as per ISO 9001:2000, ISO 17025:2005, and ISO Guide 34:2000 guidelines. Vendor identifications are proprietary, however sources of all materials used in the preparation and testing of SPEX CertiPrep CRMs are tracked and documented. For further assistance, please contact the Sales Support Department at [crmsales@spexcsp.com](mailto:crmsales@spexcsp.com).

### Instructions for Use:

Primary usage of this CRM is in neat form or diluted serially with matrix of a purity at or greater than the purity of the original matrix solution. If dilution is required the diluent must be compatible with all certified analytes and contain stabilizers appropriate for the period of intended use. The CRM can also be used as a spike or with a spike, again with appropriate compatibility considerations. All solutions should be thoroughly mixed, by shaking, prior to use and never pipetted directly from the bottle. All surfaces that come in contact with the solution must be thoroughly cleaned and leached prior to use. Dilutions should be performed only with Class A volumetric glassware.

### Method of Preparation:

Clean laboratory procedures and techniques have been used throughout the preparation. All materials, equipment, analytical instrumentation and personnel have been qualified prior to use. The highest purity acids applicable, 18 megohm, double deionized water, acid-leached triple-rinsed bottles (where appropriate), and Class A/calibrated volumetrics have been used in all preparations.

### Homogeneity:

The homogeneity of the CRM has been confirmed by procedures consistent with ISO 17025:2005, ISO Guide 34:2000, and ASTM D6362-98 Appendix X2. Random, replicate samples of the final, packaged material have been analyzed to prove homogeneity in accordance with our internal procedure 4600-HOMOGEN-1A. This is consistent with the intended use of the CRM.

### Statistical Estimator and Confidence Limits:

The certified value 'X' listed on the reverse of this document is at the 95% level of confidence and can be expressed as:

- $X = \bar{x} \pm U$  where  $\bar{x}$  = measured value,  $U$  = expanded uncertainty
  - $U = k u_c$  where  $k=2$  is the coverage factor at the 95% confidence level
- $U_c$  is obtained by combining the individual element standard uncertainty components  $u_i$ , and  $u_c = \sqrt{\sum u_i^2}$

### Certification Traveler Report:

All certified values reported were derived from the Traveler Report (SPEX CertiPrep's traceability documentation) identified by the lot number of this CRM. For further assistance, please contact the Sales Support Department at [crmsales@spexcsp.com](mailto:crmsales@spexcsp.com).

### Legal Notice:

SPEX CertiPrep reference materials are not for any cosmetic, drug or household application and are to be used only by qualified individuals who are trained in appropriate procedures. No claims against SPEX CertiPrep, Inc. of any kind whatsoever, whether based on breach of warranty, alleged negligence, or otherwise, with respect to this Reference Material shall be greater than the purchase price. In no event shall SPEX CertiPrep, Inc. be liable for any loss of profits or any incidental, special, or consequential damages.

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